

# Triple Rinse Procedure

Contact Us



**1** Fill bottle with water to 10% of its capacity.

**2** Replace cap, then shake and swirl the bottle.

**3** Pour rinse water into a labeled waste container.

**4** Repeat rinse process two more times.

**5** Deface original label and add Triple Rinse label.

**6** Place bottle in trash, recycle bin, or designated collection area.

Do not triple rinse containers with sludge, solidified material, more than residual chemicals or septum caps. Submit as chemical waste instead.

For more information regarding Lab Waste Procedures



To order Triple Rinsed lables

